



更高效 更安全 更稳定

eBlot™ L2 快速湿转仪

全新升级的湿式蛋白转印!



扫码免费试用eBlot™ L2

eBlot™ L2 是一款全新升级的蛋白快速湿转仪，能够在 5-15 分钟内同时完成 4 块小型 (mini) 凝胶或 2 块中型 (midi) 凝胶的蛋白转印过程。采用金斯瑞自主研发的技术，基于湿转的原理，集合了传统湿转稳定、均一和半干转快速、高效的优势；配套试剂不含醇，让您在清洁安全的实验环境中，高效便捷的完成转膜操作，获得一致稳定的转膜结果。



快速湿转

最快**5分钟**完成
蛋白转印



高通量

同时完成**4 mini**
或 **2 midi** 凝胶转印



安全环保

升级**试剂不含醇**
纯水稀释




适用性广

确保**一致性和可重现性**，
大、小蛋白转印

eBlot™ L2 转膜操作

01

打开转膜槽，依次平铺：
干转膜垫片-膜-凝胶-干转膜垫片

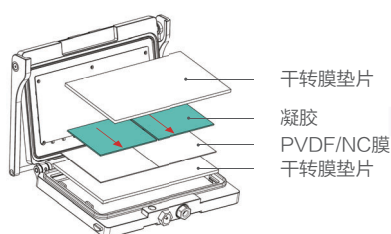

02

合上转膜槽，放在仪器对应通道中，
点击开始转膜




03

转膜结束，
抗体孵育、显影



应用数据

分别使用 eBlot™ L2 快速湿转仪和传统湿转方法进行转膜，平行对比不同分子量大小蛋白的转膜结果，从下图可以看出，eBlot™ L2 结果优于传统湿转，或者和传统湿转相当。

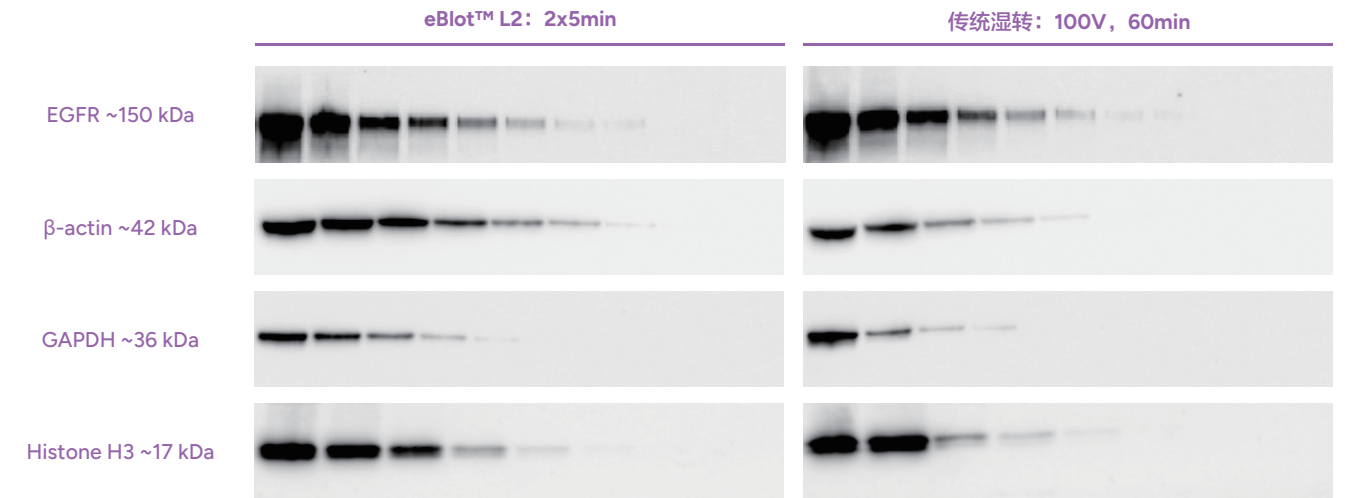

样本：HEL293 细胞裂解液；膜：NC膜，0.22 μm，凝胶：YoungPAGE™，Bis-Tris，10×8，4-12%，11 wells (GenScript, M00929)
上样：Lane 1~8: 样品 20ug, 10ug, 5ug, 2.5ug, 1.25ug, 0.62ug, 0.31ug, 0.15ug (EGFR); Lane 1~9: 样品 40ug, 20ug, 10ug, 5ug, 2.5ug, 1.25ug, 0.62ug, 0.31ug, 0.15ug (β-actin, GAPDH, Histone H3)
一抗：Recombinant Anti-EGFR antibody, mAb, rabbit (abcam, ab52894), 1:1000; THE™ beta Actin antibody, mAb, mouse (GenScript, A00702), 1:5000; GAPDH antibody, mAb, mouse (GenScript, customized), 1:1000; Histone H3 antibody, pAb, rabbit (GenScript, customized), 1:1000
二抗：Anti-mouse IgG (HRP) (GOAT) antibody peroxidase (Rockland, 610-103-121), 1:10000; Anti-rabbit IgG (HRP) (GOAT) antibody peroxidase (Rockland, 610-103-122), 1:10000

产品列表

产品名称	规格	产品货号
eBlot™ L2 快速湿转仪	1 unit	L00981C
eBlot L2 转膜浓缩试剂盒, midi	1 kit	L01015C-75
SurePAGE™ Midi, Bis-Tris, 4-12%, 20/26 wells	10 PK/BOX	M00995 / M00996
SurePAGE™ Midi, Bis-Tris, 12%, 20/26 wells	10 PK/BOX	M01110 / M01111
SurePAGE™ Midi, Bis-Tris, 10%, 20/26 wells	10 PK/BOX	M01106 / M01107

2hrs Western Blot 流程

