

Green *Taq* DNA Polymerase

Cat. No.: E00043

Overview

Description	Green Taq DNA Polymerase is designed to increase the stability of the Taq enzyme for more convenient transport and applications. The polymerase (1000 U) is designed for 400 rxns if 2.5 U are used per 50 ul PCR reaction. It can be stable for six months if stored at 4°C or for one month without significant loss of activity at ambient temperature. Our technology can increase the stability of Green Taq DNA at higher temperature (72°C). The higher yield PCR product can be harvested in long PCR amplification.
Key Features	High Stability: The enzyme remains stable for more than six months when stored at 4°C or for one month when stored at room temperature (25°C) without significant activity loss. High PCR Yield: The Green Taq DNA polymerase has longer enzyme half-life and therefore increases the PCR yield when amplifying long DNA. Terminal Transferase Activity: Taq DNA polymerase has terminal transferase activity that results in the addition of a single nucleotide (adenosine) at 3' -end of the extension product. Stop refrigerating and save energy.

Properties

Unit Definition	One unit is defined as the amount of enzyme that can incorporate 10 nmol of dNTP into acid-insoluble material in 30 minutes at 74°C.
Formulation	The polymerase is designed for 400 rxns if 2.5 U are used per 50 ul PCR reaction. It can be stable for six months if stored at 4°C or for one month without significant loss of activity at ambient temperature.
Reaction Buffer	PCR Reaction Buffer (with Mg ²⁺): 100 mM Tris-HCl, 15 mM MgCl ₂ , 500 mM KCl and 1% Triton X-100, pH 8.3 (4°C).
Storage & Stability	This product is shipped at ambient temperatures for several weeks. Green Taq DNA Polymerase will remain stable for six months if stored at 4°C and for two years if stored at -20°C. Please refer to the product label for the specific shelf life of each batch.

Applications

The applications of Green Taq DNA Polymerase include the following:

- PCR* (For simple templates, this enzyme can be optimized for amplification of PCR products up to 10 kb; However, for complex templates, this enzyme can be used for amplification of PCR products up to 3 kb.)
- 3' A-tailing of blunt ends (T/A-cloning)
- Primer extension
- DNA labeling reactions

For research use only. Not intended for human or animal clinical trials, therapeutic or diagnostic use.

Manufacturer: Nanjing GenScript Biotech Co., Ltd. No. 28Yongxi Road, Jiangning District, Nanjing, Jiangsu, China